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Instructions

TPP tissue culture flasks are intended for cell culture with necessary sterile handling techniques. The growth area is opto-mechanically treated for optimal cell adhesion and proliferation. The angled neck decreases the risk of media contact with the inside of the screw cap.

- Open the product packaging and inspect for intact bag and product.
- Fill medium: Introduce medium or inoculate cells with a pipette alongside the wall, or direct injection on the growth area of the flask.
- Tightly screw closed the filter screw cap with the integrated hydrophobic membrane for sterile gas exchange
- Incubate: Lay the flask on the incubator shelves for short or long term storage.
- **Open re-closable flask:** Complete access from above to flask chamber with multi-use re-closable liquid-tight lid.
- Harvest cells with large access from above to flask chamber
 - Harvest adherent cells by trypsinization or scrape with the cell scraper.
 - Harvest suspension cells by exchange of medium.

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Advice and Recommendation

*Stable ring enables slip-free stacking of multiple flasks also in combination with the standard TPP tissue culture flasks.
 Barrier creates a growth area of 115 cm².
 Both flasks are available with Filter screw cap with a hydrophobic PTFE membrane with a pore size of 0.22 µm only. If a «VENT» screw cap is used, the build-up of inside pressure of more than 0.03 bar could press open the flask.
 A controlled level base surface and regular use of feeding medium are essential to uniform tissue growth.
 Sources of vibrations must be eliminated to avoid circular growth pattern.*

Technical Data

Materials

Flask	PS
Screw cap	PE
Membrane	PTFE, pore size 0.22 µm
Gasket	non-toxic gasket

Abmessungen	Measurement	90552	90652
Wiederverschliessbar	<i>Re-closable</i>	✓	✓
Steg	<i>Barrier</i>	---	✓
Länge mm	<i>Length mm</i>	210	210
Breite mm	<i>Width mm</i>	122	122
Höhe mm	<i>Heights mm</i>	50	50
Wachstumsfläche cm ²	<i>Growth area cm²</i>	115	115
empf Volumen ml	<i>Recom. volume ml</i>	15 – 45	15 – 33
Öffnung mm	<i>Opening mm</i>	105 x 105	105 x 105
max Volumen ml	<i>Max volume ml</i>	100	100